

Emergency Eyewash Station Weekly Inspection Checklist

How to Test:

- Is the area around the station clear of obstructions?
- Are the eyewashes clean and free of dirt/debris?
- Flush eyewash for 30 seconds weekly
- Is the water clear?
- Is the plumbing free from leakage?
- Do the covers/caps open without assistance?
- Are the jets working properly
(both spigots are delivering evenly, and with sufficient but not too great of a flow)?

How to obtain assistance:

Questions, maintenance requests, training information
Environmental Health and Safety: 4-5084 or
ehs_labsafety@utk.edu

Maintenance requests may also be called in directly
Facilities Services (Zone Maintenance): 4-5346

Emergencies:

Hold eyes open and flush for at least 15 minutes. CALL 911
and seek medical assistance

Date	Initials	Comment	Date	Initials	Comment	Date	Initials	Comment