

Emergency Eyewash Station Weekly Inspection Checklist

How to Inspect & Test:

- Is the area around the station clear of obstructions?
- Is the eyewash clean and free of dirt/debris?
- Flush eyewash for 30 seconds weekly
- Do the covers/caps open without assistance?
- Is the plumbing free from leakage?
- Are the jets working properly (both spigots are delivering evenly, and with sufficient but not too great of a flow)?
- Is the water clean and clear?

How to obtain assistance:

Questions, maintenance, training
 EHS: 865-974-5084 | ehs_labsafety@utk.edu

Maintenance
 requests may also be called in directly
Facilities Services: 865-946-7777

Emergencies:
 Hold eyes open and flush for at least 15 minutes.
 CALL 911 and seek medical assistance

Date	Initials	Comment		Date	Initials	Comment		Date	Initials	Comment